

## Splitting hPSCs with dispase


Mark Tomishima, Sloan-Kettering Institute, 430 East 67th Street New York, NY 10065 US

### Introduction

This protocol outlines the techniques used to routinely passage hPSCs in the SKI Stem Cell Research Facility at Sloan-Kettering. We prefer to culture our cells on mouse embryo fibroblasts (MEFs) but occasionally grow the cells feeder free for particular applications, such as nucleofection, viral transduction or karyotyping.

We use Dispase for all applications requiring colony passage including feeder-dependent and independent expansion as well as for cryopreservation.

### Flow Chart

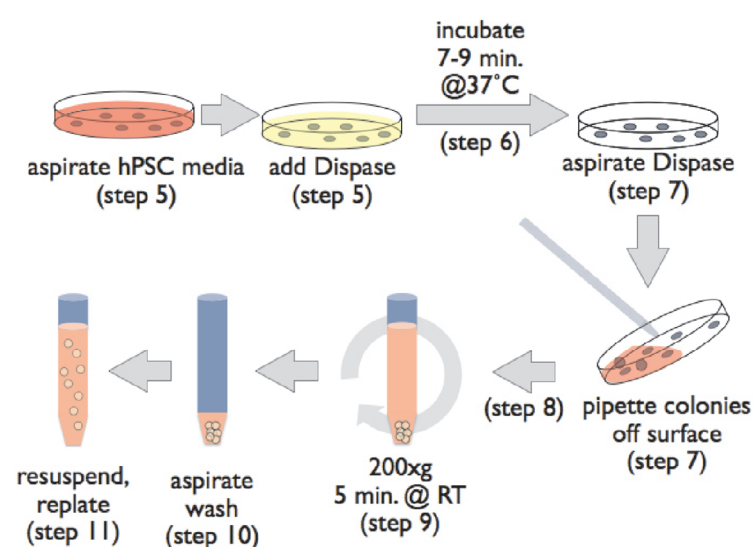

### Matrigel coating dishes

1. To expand cells feeder free, you must first prepare Matrigel-coated dishes. The preferred method for doing this is to remove a 1 ml aliquot (in a 50 ml conical stored at  $-20^{\circ}\text{C}$ , see Materials) and place on ice overnight. In a pinch, you can warm it carefully in your hand, but make sure to place it on ice before it completely melts.
2. Add 19 ml of cold DMEM:F12 to the Matrigel. Sometimes, small pieces of polymerized Matrigel form. These do not interfere, but you can filter them out using a  $100\ \mu\text{M}$  strainer if desired (see Materials).

Last revised March 28, 2012. Published June 10, 2012. This chapter should be cited as: Tomishima, M., Splitting hPSCs with dispase (June 10, 2012), StemBook, ed. The Stem Cell Research Community, StemBook, doi/10.3824/stembook.1.80.1, <http://www.stembook.org>.

**Copyright:** © 2012 Mark Tomishima. This is an open-access article distributed under the terms of the Creative Commons Attribution License, which permits unrestricted use, distribution, and reproduction in any medium, provided the original work is properly cited.

**Correspondence:** Mark Tomishima. E-mail: [tomishim@mskcc.org](mailto:tomishim@mskcc.org)

### Splitting hPSCs with dispase

3. Add Matrigel to your surface. Examples of volumes by surface area/vessel type:
4. Incubate in the hood for one hour.

#### Passage of cells using Dispase

5. When cells are ready for passage, remove hPSC media and add the appropriate volume of 5 mg/ml Dispase.
6. Incubate for 7–9 minutes at 37°C in the incubator. After 7 minutes, check on the microscope to verify that the edges have started to detach. If they have not, return to the incubator for another minute and then check again.
7. Once they have started to peel, remove the Dispase and detach colonies from the cell surface with a small volume of hPSC media. Generally, this can be the same volume that you used for the Dispase (e.g., 2 ml for a 6 cm dish). You should be able to dislodge the colonies with relatively gentle pipetting. Try to keep the colonies as large as possible while removing them.
8. Add the detached colonies to an adequate volume of hPSC media, DMEM:F12 or even PBS to wash. Dispase is not inhibited with media, so dilution is required to inactivate the enzyme. I prefer to use 5 times the volume of dispase used per wash.
9. Spin at 200×g for 5 minutes at room temperature. Alternatively, you can just let the clusters settle with gravity for 5 minutes in the hood. This might help to eliminate single cells that are thought to contribute to karyotype problems in hPSC cultures.
10. Aspirate wash, carefully avoiding cell clusters. (Wash a second time if you are plating without feeders)
11. Resuspend in hPSC media. I typically resuspend in 12 ml of hESC media and plate 1 ml to a new dish (1:12), but this must be adjusted for each culture. Generally it is in the range of 1:6–1:20.

#### Troubleshooting

Use the Dispase specified in Materials: not all Dispases are the same. This protocol will only work for this preparation. We have used Worthington in the past too, but it is stronger and requires a second wash. When passaging cells without feeders, they are exquisitely sensitive to any residual enzyme left with the cells. Make sure you wash with a large volume and wash 2 times. If you are having trouble with the cells not attaching, you might even want to wash 3 times.

Keep in mind that as long as the colonies are still attached, you can wash the colonies by adding media/PBS to the dish and aspirating in place. This is much faster than spinning them down. At least one wash/spin should be done after you have detached colonies, however.

Dispase can not be inhibited by media so it is not necessary to use complex media to wash cells. Dilution is the only way to eliminate the enzyme.

#### Preplating to remove residual feeders

If you are moving your hPSCs from feeders to a feeder independent condition, you might want to preplate cells to remove feeders. Basically, you prepare your cells as you normally would be let them attach to gelatin-coated tissue culture treated plastic for one hour in the incubator. Under these conditions, the MEFs will adhere to the dish but the hPSCs will poorly attach to this plastic. After the hour incubation, remove the dish, pipette over the surface of the dish gently to dislodge the hPSCs, and transfer to a new Matrigel-coated dish leaving the attached fibroblasts behind.

#### Materials

Reagents	Supplier	Catalog#	Notes
Dispase 5 mg/ml	StemCell Technologies	07913	Do not substitute!
Matrigel	BD Bioscience	352340	We only use lots over 10 mg/ml protein
45 µm nylon mesh strainer	BD Falcon	352340	
FGF2	R&D Systems	233-FB-001 MG/CF	Dissolve in PBS with 0.1% BSA to 100 ug/ml

---

Splitting hPSCs with dispase

---

hPSC media

800 ml DMEM/F12	Life Technologies	11330-032
200 ml KSR	Life Technologies	10828-028
5 ml L-glutamine	Life Technologies	25030-081
10 ml MEM NEAA	Life Technologies	11140-050
1 ml 2-mercaptoethanol	Life Technologies	21985-023

Sterile filter, then add FGF2 to 6 ng/ml. Use within 2 weeks.